Bestell-Nr. ZS - 025
Order no. ZS - 025

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Function

Input:
-PTC sensor
-open door alarm, normally closed (optional)
-shaker failur alarm, normally closed (optional)

Output:
- collective failur indicator, external alarm
- temperature 1 alarm
- temperature 2 alarm
- disconnection of relais cooling system when low temperaturalarm goes off

Serial Interface RS 485 (optional)

Alarmdisplay:
- door open (optional)
- shaker off (optional)
- alarm temp 1
- alarm temp 2 (option reversierbar),
- accu failure
- mains off

Mains off
- lighting of the alarm display is blinking.
- display indicated mains off
- cooling regulator off, no display .

Standard adjustments:
- overtemp T1, undertemp T2 (cooling off)

Option
- overtemp T1, overtemp T2

Battery (accu) usage:
- approximately 4 hrs. Alarmdisplay during a power outage, when accu fully charged.
- charging time 48 hrs.

Button Functions

SET
When pressing the SET button, the alarm sensor present value of the actual temperature will be displayed.
TEST
When pressing the TEST button, a diagnostic program of the alarm panel will be started. as long as the TEST button is pressed, a hooter will sound.

Display: alarm control (Software number)
When pressing QUIT and SET at the same time during the test, the program level will be reached. During an alarm display, the program level can be reached by pressing the QUIT and SET buttons simultaneously. After the display "alarm off", release first QUIT, then SET. Now the program data can be changed.

Arrow Keys Change values by using the arrow keys

Alarm

Alarms are shown on the display in a numerical manner and are signalised acoustically by a hooter.

Display of the alarm after at least 30 seconds.
Collective failure indication relay is connected delayed. Can be set individually.
During an alarm, the relay is shut off. In case of an alarm, the hooter will sound simultaneously.

It can be receipted by pressing the QUIT/TEST button. Even if the alarm was not receipted (visual display), the hooter will sound during a new alarm.

The QUIT/TEST button must be pressed longer than **20 seconds to delete** the alarm display when the alarm has ended.
The alarm display cannot be receipted if there is another alarm to come. By using the QUIT/TEST button, a receipt over the failure indication to the external alarm display is printed.

Door open
- Acoustic signal in intervals after 3 minutes (adjustable). **No alarm exit.**
- Display: door open
- **Alarm can only be deleted by pressing the QUIT button (approx. 20 sec.)**
- Delayed alarm-output (after adjusted collective alarm delay-time)
- After elapsed time: Display: door open (blinking), permanent sounding of hooter.
- **Alarm can only be deleted by pressing the QUIT button (approx. 20 sec.)**

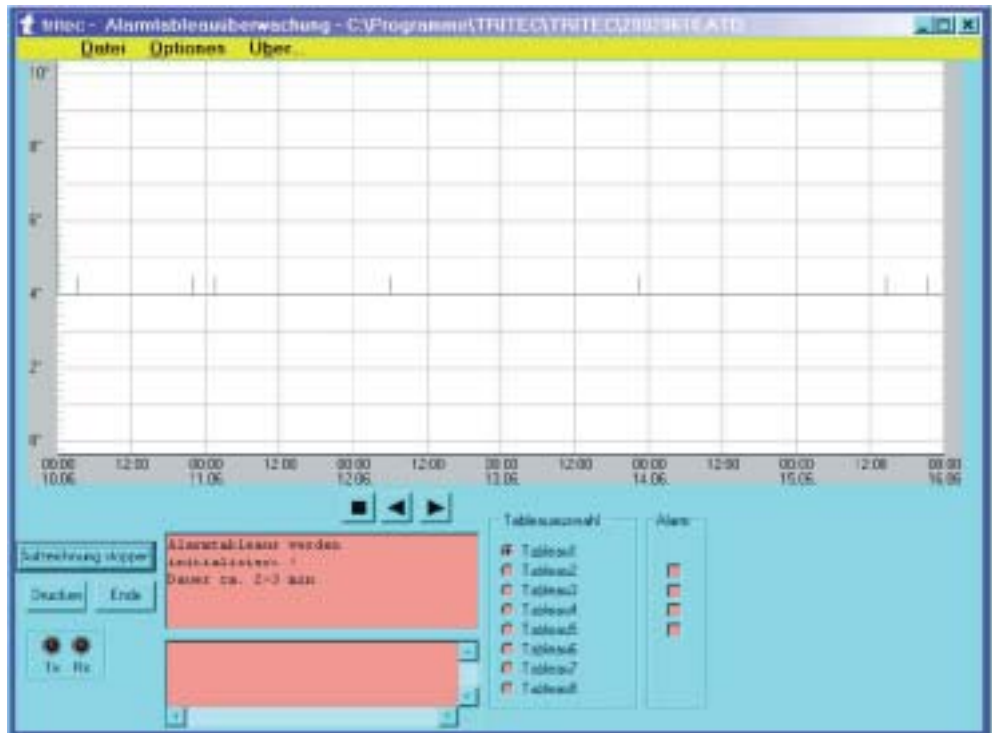
Shaker off (optional)
- As soon as the contact between the shaker and the alarm panel board is cut off, an alarm will sound
- delayed collective alarm-relay (adjustable)
- **Alarm can only be deleted by pressing the QUIT/TEST button (approx. 20 sec.)**

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Monitor description of graphic records (800 x 600 pixel).
Display of the connected panels (cabinets), fault signalling display, adjustable scaling, data files weekly.

Description:
- weekly Display (raster 12 hours)
- increase the raster to 1 hour,
- increase the raster to 5 minutes

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Printout of the records in the shown zoom-area (DinA4-landscape format). The printout also includes the appropriate alarm journal.

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